

# User Manual

Kohler Illumination  
Binocular Compound  
Microscope

Model M828D5



[MicroscopeNet.com](http://MicroscopeNet.com)

# Table of Contents

i. Caution.....	1
ii. Care and Maintenance.....	2
1. Component Illustration.....	3
2. Installation.....	4
3. Operation.....	6
4. Specifications.....	8
5. Troubleshooting Guide.....	9

## i. Caution

1. Find the “UP” sign and place the Styrofoam container on your table or bench so that the arrow is pointing upward. Open the shipping carton carefully to prevent any accessory items (i.e. objectives or eyepieces) from dropping and being damaged.
2. Do not discard the molded Styrofoam container. The container should be retained should the microscope ever requires reshipment.
3. Keep the instrument out of direct sunlight, high temperature or humidity, and dusty environments. Ensure that the microscope is located on a smooth, level and firm surface.
4. If any specimen solutions or other liquids splash onto the stage, objective or any other component, disconnect the power cord immediately and wipe up the spillage. Otherwise, the instrument may be damaged.
5. **Important:** the lamp, lamp housing and adjacent parts will become very hot. Do not touch these parts until they have completely cooled. Never attempt to handle a hot halogen bulb.
6. All electrical connectors (power cord) should be inserted into an electrical surge suppressor to prevent damage due to voltage fluctuations.
7. For safety when replacing the halogen lamp or fuse, be sure the main switch is off, unplug the power cord, and only replace the halogen bulb after the bulb and the lamp house has completely cooled.
8. Confirm that the input voltage indicated on your microscope corresponds to your line voltage. The use of a different input voltage other than that as indicated will cause severe damage to the microscope.

## ii. Care and Maintenance

1. Do not attempt to disassemble any component including eyepieces, objectives or focusing assembly.
2. Keep the instrument clean; remove dirt and debris regularly. Accumulated dirt on metal surfaces should be cleaned with a damp cloth. More persistent dirt should be removed using a mild soap solution. **Do not use organic solvents for cleansing.**
3. The outer surface of the optics should be inspected and cleaned periodically using an air stream from an air bulb. If dirt remains on the optical surface, use a soft cloth or cotton swab dampened with a lens cleaning solution (available at camera stores). All optical lenses should be swabbed using a circular motion. A small amount of absorbent cotton wound on the end of a tapered stick makes a useful tool for cleaning recessed optical surfaces. Avoid using an excessive amount of solvents as this may cause problems with optical coatings or cemented optics or the flowing solvent may pick up grease making cleaning more difficult. Oil immersion objectives should be cleaned immediately after use by removing the oil with lens tissue or a clean, soft cloth.
4. Store the instrument in a cool, dry environment. Cover the microscope with the dust cover when not in use.



## 2 Installation

### 2.1 Installation of the binocular viewing head

- 1) Loosen the thumb screw (11) on the top of the body (12) and remove the plastic cover on the top.
- 2) Remove the cap on the circular dovetail of the binocular viewing head (10)
- 3) Insert the dovetail into the socket on the top of body (12); ensure that it is completely seated into the socket; tighten the thumb screw (11).



**Caution:** Do not release the head from your hand grip until you are sure the head is installed securely.

### 2.2 Installation of the eyepieces

- 1) Remove the protective caps from the eyepiece tubes (3).
- 2) Insert the eyepieces (1) into the eyepiece tubes (3).

### 2.3 Installation of the objectives

- 1) Adjust the coarse focus knob (14) until the mechanical stage (7) is at its lowest position.
- 2) Install the 4x objective into the nosepiece (4). Then in a clock-wise direction, rotate the nosepiece and install each succeeding higher magnification objective.

**Note:** Inspect the objectives frequently for dirt or oil; clean if necessary. Use the 10x objective to initially focus the image of your specimen. When changing the objective magnification, rotate the objective nosepiece until you hear a “click” sound. This ensures the objective is centered in the optical light path.

### 2.4 Installation of the filter

- 1) Swing out the filter holder under the condenser (21).
- 2) Put the filter onto the holder, swing the holder back in.



### 2.5 Installing (or changing) the halogen bulb

- 1) Turn the power off and disconnect the power cord.
- 2) Allow some time to cool down the lamp.
- 3) Turn over the microscope on its side; find the bulb compartment at the bottom.
- 4) Open the cover of the bulb compartment by loosening the thumb screw. Take out the dead bulb and insert the new bulb. Be sure the pins on the bulb are completely inserted into the lamp socket. You may also loosen the two screws on the cover to adjust the position of the bulb to get centered and even brightness. Screw the cover on.

**Caution:** Before you turn over the microscope, be sure to take the eyepieces off and be certain that the head is securely locked by the thumb screw.



### 2.6 Replacing the fuse

- 1) Turn off the power and disconnect the power cord.
- 2) Turn over the microscope on its side; find the fuse holder at the bottom of the base.
- 3) Turn the fuse holder counter-clockwise to take it out, replace the fuse, and then turn

it on clockwise.

**Caution:** Before you turn over the microscope, be sure to take the eyepieces off and be certain that the head is securely locked by the thumb screw.

### 2.7 Installing the mirror (optional, your model may not have one)

- 1) Unplug the power cord.
- 2) Thread off the field diaphragm on the base.
- 3) Thread the black disc onto the base and then insert the mirror into the hole at the center of the black disc. You may try to get reflected ambient light on either side of the mirror with different angles for best result.



**Note:** The mirror is only used when there is a power outage or you are in the field and no power available.

## 3 Operation

### 3.1 Adjusting illumination

- 1) Connect the power cord and turn on the main switch (18).
- 2) Turn the variable intensity dial (17) to increase or decrease the brightness.

### 3.2 Placing specimen

- 1) Place the slide on the mechanical stage (7).
- 2) Use the slide holder (6) to gently secure the slide.
- 3) Turn the X and Y translational knobs (16) to position the specimen for viewing.

**Caution:** Be sure not to allow an objective to touch a specimen slide when changing objectives.

### 3.3 Adjusting interpupillary distance

While observing with both eyes, hold the left and right eyepiece tubes (3) then slide the tubes in and out. The interpupillary distance is correct when the left and right fields of view converge completely into one image.

### 3.4 Adjusting eyepiece diopter

- 1) Rotate the 10x objective into position.
- 2) Rotate the diopter ring (2) on the right eyepiece tube (3) until its numerical value is the same as your interpupillary distance, for example, 65 in the right figure.
- 3) Close your left eye and bring the specimen into focus following the focusing procedures in 3.5.
- 4) Close your right eye and bring the same specimen into clear sharp focus by adjusting the diopter ring (2) on left eyepiece tube (3) only. Do not use focus knobs at this step.
- 5) Since both sides are adjustable, you may also do the above in the opposite way, in other words, left eye first and right eye second.



### 3.5 Focusing

- 1) With the 10x objective in position, raise the mechanical stage (7) using the coarse focus knob (14) until the specimen is close to the objective
- 2) Turn the coarse focus knob (14) until the specimen is in focus.
- 3) Use the fine focus knob (15) to obtain a sharp image.
- 4) You may now switch to another magnification objective.

**Tips:** To prevent your specimen slide from making contact with an objective, raise the stage to its highest position without contacting the 100x objective, then tighten the focus stop lever.

### 3.6 Adjusting condenser

- 1) Turn the condenser focus knob (20) to raise or lower the condenser (22).
- 2) Raise the condenser (22) when using high power objectives and lower it when using low power objectives.

**Note:**

- The centering of the condenser and the light axes of the objective are factory adjusted. Do not attempt re-adjustment.
- The highest position of the condenser has been factory adjusted. Do not attempt re-adjustment.

### **3.7 Adjusting Kohler illumination**

- 1) Connect the power cord and turn on the main switch.
- 2) Have the specimen in focus.
- 3) Close the field diaphragm (24) so that the edges of the field diaphragm show in the field of view.
- 4) Turn the condenser focus knob (20) to make the image of field diaphragm edges clear
- 5) Adjust the centering screws (23) to make the image of field diaphragm edges in the view center.
- 6) Open the field diaphragm (24) so that its edges are just out of the view field.
- 7) Adjust the aperture iris diaphragm lever (19) to get proper contrast of the specimen.
- 8) Adjust the intensity dial (17) to get proper light intensity.

### **3.8 Adjusting focus tension**

The tension has been pre-set at the factory. If the mechanical stage drops by itself, rotate the tension adjustment collar (13) located inside the focus knob (14) on the power switch side until the tension is maintained.

## 4 Specifications

### General

Model	M828D5
Total Magnification	40X, 64X, 100X, 160X, 200X, 320X, 400X, 640X, 1000X, 1600X
Viewing Head	Binocular, inclined 45°, swiveling 360°
Interpupillary Distance	Sliding adjustment, 55mm ~ 75mm
Diopter Adjustment	On both eyepiece tubes
Eyepieces	1 pair of WF10X 1 pair of P16X
Nosepiece	Revolving, quintuple
Objectives	Plan objectives 4X, 10X, 20X, 40X(spring), 100X(spring, oil)
Condenser	Abbe, NA=1.25, w/ iris diaphragm and filter holder Rack and pinion adjustment
Focus Mechanism	Coaxial coarse and fine focusing knobs on both sides w/ focus stop
Stage	Double layer mechanical stage Dimension: 140mmx140mm (5-1/2" x 5-1/2") Movement range: 75mm X 50mm (3" x 2")
Illumination	Transmitted Kohler illumination system, adjustable center of field diaphragm Light bulb: 6V/20W, Halogen, Variable intensity
Power Supply	AC 100V ~ 240V, 50/60HZ (US and Canada plug)
Dimension	28cm x 19cm x 40 cm (11" x 7-1/2" x 15-3/4")
Net weight	5.34kg (11lb 12oz)
Package weight	6.6 kg

### Magnifications

Eyepiece	10X	16X	10X	16X	10X	16X	10X	16X	10X	16X
Objective	4X		10X		20X		40X		100	
Magnification	40X	64X	100X	160X	200X	320X	400X	640X	1000X	1600X

## 5 Troubleshooting Guide

### OPTICAL PROBLEMS

<b>Problem</b>	<b>Cause</b>	<b>Solution</b>
Darkness at the periphery or uneven brightness in the field of view	Revolving nosepiece not in click stop position	Revolve the nosepiece to click-stop position by swinging the objective correctly into the optical path
	The light source of the bulb is not at the center	Adjust the position of the bulb
Dirt or dust on the view	Dirt or dust on the lens eyepiece, condenser, objective, collector lens or specimen	Clean the lens with a camera cleaning kit
	Dirt or dust on the specimen	Clean the specimen
Poor image quality	No slide cover attached to the slide	Attach a 0.17mm slide cover
	Slide cover is too thick or thin	Use a slide cover of the appropriate thickness (0.17mm)
	Slide may be upside down	Turn slide over so the cover-glass faces up
	Immersion oil is on a dry objective (especially the 40x)	Check the objectives, clean if necessary
	No immersion oil used with 100x objective	Use immersion oil
	Air bubbles in immersion oil	Remove bubbles
	Condenser aperture is closed or open too much	Open or close properly
	Condenser is positioned too low	Position the condenser upward
	Kohler illumination system not set properly	Set up the Kohler illumination system according 3.4 on page 8.

### ELECTRICAL PROBLEMS

<b>Problem</b>	<b>Cause</b>	<b>Solution</b>
Lamp does not light when switched on	No electrical power	Check power outlet Check power cord connection
	Lamp bulb burnt out	Replace bulb
	Fuse blown out	Replace fuse

**IMAGE PROBLEMS**

<b>Problem</b>	<b>Cause</b>	<b>Solution</b>
Image moves while focusing	Specimen rises from stage surface	Secure the specimen in the slide holder
	Revolving nosepiece is not in the click-stop position	Revolve the nosepiece to the click-stop position
Image tinged yellow	Blue filter not used	Use daylight blue filter
	Lamp intensity is too low	Adjust the light intensity by rotating the intensity control dial
Image is too bright	Lamp intensity is too high	Adjust the light intensity by rotating the intensity control dial
Insufficient brightness	Lamp intensity is too low	Adjust the light intensity by rotating the intensity control dial
	Aperture diaphragm closed too far	Open to the proper setting
	Condenser position too low	Position the condenser upward

**MECHANICAL PROBLEMS**

<b>Problem</b>	<b>Cause</b>	<b>Solution</b>
Image will not focus with high power objectives	Slide upside down	Turn the slide over so the cover glass faces up
	Cover glass is too thick	Use a 0.17mm cover glass
High power objective contacts slide when changed from low power objective	Slide upside down	Turn the slide over so the cover glass faces up
	Cover glass is too thick	Use a 0.17mm cover glass
	Diopter adjustment is not set properly	Readjust the diopter settings
Slippage of focus when using the coarse focusing knob. Fine focus is ineffective	Tension adjustment is set too low	Increase the tension on the focusing knobs
	Tension adjustment is set too high	Loosen the tension on the focusing knobs